



ImmEdge™ PEN

Catalog Number H-4000

Description: Hydrophobic barrier pen for immunohistochemistry and *in situ* hybridization. ImmEdge™ Pen provides a heat-stable water-repellent barrier that keeps reagents localized on tissue sections.

Unit Size: 2 pen set

Storage Conditions: Room temperature with pen tip down and tightly capped.

INSTRUCTIONS FOR USE

1. Shake vigorously with lid on.
2. Depress tip to saturate, if necessary.
3. Circumscribe tissue section with the ImmEdge™ Pen.
4. Replace pen cap. Store pen tip down.
5. Allow ImmEdge™ Pen reagent to dry before immersing slides in aqueous solutions.

DISPOSAL OF REAGENTS

Dispose in an approved landfill.

MSDS available online at: www.vectorlabs.com

For Professional Use Only.

Revised 01/18