



VECTOR®
NUCLEAR FAST RED
COUNTERSTAIN
Catalog Number H-3403

This nuclear counterstain is designed to be used after completion of immuno-histochemical staining or for routine histology. Nuclei in stained sections will be pink to light red.

INSTRUCTIONS FOR USE:

1. Rinse slides in tap water.
2. Immerse slides in Vector® Nuclear Fast Red counterstain or apply counterstain directly to slide, completely covering the tissue section.
3. Incubate sections for 1 to 10 minutes to obtain desired stain intensity.
4. Wash slides in tap water for 10 minutes. **Washing for longer than 10 minutes may lead to decreased counterstain intensity.**
5. Dehydrate, clear, and mount.

NOTES:

Counterstains should be optimized for each tissue type, antigen unmasking protocol, and immunocytochemical staining intensity desired.

A precipitate may form in this product over time. This will not affect staining efficacy. If preferred, product can be filtered immediately before use.

See reverse side for substrate compatibility.

Use ordinary precautions to avoid contact with skin and eyes.

Store at room temperature.

This product is for research use only.