


VECTASHIELD® Antifade Mounting Medium

Cat. No.: H-1000

Storage: Refrigerate

Refractive Index: 1.45

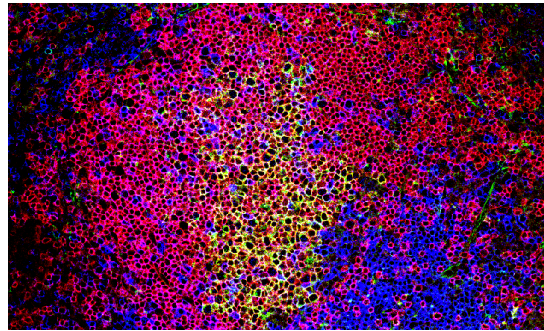

Germinal center (GC) after acute viral infection. T cells (in blue) B cells (in purple), GCs (in green). Mounted in VECTASHIELD Antifade Mounting Medium.

DESCRIPTION

VECTASHIELD Antifade Mounting Medium is a unique, stable formula for preserving fluorescence. VECTASHIELD Antifade Mounting Medium prevents rapid photobleaching of fluorescent proteins and fluorescent dyes (see website for compatibility).

Each lot is tested for photobleaching retardance with qualitative and quantitative methods.

APPLICATION

To mount tissues or cells on a slide dispense VECTASHIELD Mounting Medium onto the specimen using a disposable pipet or pipet tip. Small drop volumes of approximately 25 ul (per 22 mm x 22 mm coverslip) are recommended. Coverslip the specimen and allow VECTASHIELD Mounting Medium to disperse over the entire section. Tightly seal the bottle with the screw cap after use and store in an upright position.

VECTASHIELD Antifade Mounting Medium does not solidify, but remains a liquid on the slide and can be stored without sealing. If desired, coverslips can be sealed around the perimeter with nail polish or a plastic sealant. Mounted slides should be stored at 4 °C, protected from light.

VECTASHIELD Antifade Mounting Media Products:

Product	Counterstain	Cat. No.	Unit Size	Hardening
VECTASHIELD Antifade Mounting Medium	none	H-1000	10 ml	no
	DAPI	H-1200	10 ml	no
	PI	H-1300	10 ml	no
VECTASHIELD HardSet™ Antifade Mounting Medium	none	H-1400	10 ml	yes
	DAPI	H-1500	10 ml	yes
	TRITC-Phalloidin	H-1600	10 ml	yes
VECTASHIELD Vibrance™ Antifade Mounting Medium	none	H-1700	2 ml, 10 ml	yes
	DAPI	H-1800	2 ml, 10 ml	yes